


TWO STAGE COMPRESSOR TEST SET F865

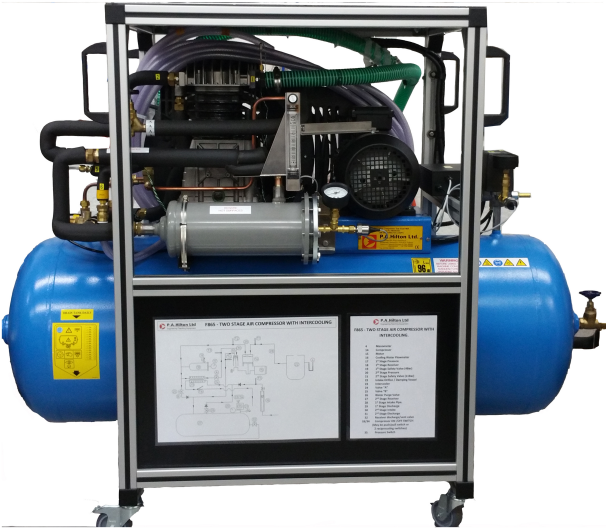

Year 1
study

3 to 4
participants

Space required
2x1.5x2m

Features

- Allows Investigation of a Single and Two Stage Compressor (with, or without intercooling) at a Range of Delivery Pressures.
- Safe and Suitable For Student Operation.
- Instrumentation Allows Detailed Analysis of Compressor Performance.
- Optional Computerised Data Acquisition Upgrade

Description

The reciprocating air compressor is a fundamental subject for thermodynamic analysis and study. The process of induction and compression is a cyclic one. However by the use of a large reservoir on intake and discharge the system can be examined as a steady flow process.

A floor standing unit with belt driven two stage reciprocating air compressor that may be operated as a

single stage, two stage, or two stage with and without intercooling system. Instrumentation allows the drive power to be measured together with the motor speed.

Compressor speed is measured through the fixed pulley ratio between the drive motor and compressor. The intake air passes through an orifice for flow measurement to a vessel that reduces intake pulsations and allows various intake conditions to be established. Valves allow an intercooler between the low and high pressure stages to be included or removed from the circuit to demonstrate and investigate comparative performance.

Valves also allow the unit to be operated in single stage or two stage mode with and without intercooling. The 1st stage and 2nd stage pressures are indicated on gauges and all of the relevant system temperatures are recorded by thermocouple sensors connected to multiple selector switches and a digital temperature indicator.

The flow measurement orifice is connected to a panel mounted manometer to allow differential pressure measurement. Air is delivered to a receiver fitted with a high pressure cut out and relief valve for operator safety. A throttle valve allows discharge pressure to be controlled and adjusted.

An optional Computerised Data Acquisition system **FC865A** is available complete with its own Installation, Operating and Maintenance Manual. The F865 has been designed so that the upgrade components can be fitted at any time during the life of the unit by any competent technician.

Alternatively the F865 can be ordered with the Computerised Data Acquisition system already fitted.

Related laws

- Thermodynamics
- Energy Conservation
- Mechanical Engineering
- Plant and Process Engineering
- Automotive Engineering
- Fluid Mechanics

Learning capabilities

- Investigation Of Variation In Air Flow Rate, With Compressor Pressure Ratio:- With And Without Intercooling
- Investigation Of Variation In Volumetric Efficiency With Compressor Pressure Ratio:- With And Without Intercooling
- Investigation Of Variation Of Isothermal Efficiency With Compressor Pressure Ratio:- With And Without Intercooling
- Investigation Of The Compressor Performance Relative To Electrical Power, Shaft Power, And Heat Loss:- With and Without Intercooling

Technical Specification

- Working Pressure: max 10bar gauge
- Temperature Range: ambient – 180°C approx.
- Motor power: 3kW
- Pressure vessel capacity: 200L

- Two Cylinder Compressor
- Low Pressure Cylinder: Bore 95mm; Stroke 50mm
- High Pressure Cylinder: Bore 50mm; Stroke 50mm
- Six temperature Measuring points
- Noise levels measured around the apparatus. (*These will vary dependent upon specific installation. Measurement parameters: 0.5m @45°): at operator's ear level 100.2db; in front of the apparatus 96.3db; back the apparatus 102.5db; left-hand side of the apparatus 97.9db; right-hand side of the apparatus 98.4db; at exhaust of the vessel. 106.3db

Recommended Ancillaries

- FC865A – Computerised Data Acquisition System

What's in the Box?

- 1 x F865
- 2 x 5m Cooling water supply and drain hose
- 1 x Manometer Fluid
- 1 x 3m Manometer hose
- 1 x Discharge pipe
- 2 x Ear Defenders
- 2 x Safety spectacles
- 2 x Pair of gloves
- Instruction manual
- Packing List
- Test Sheet
- Wiring Diagram

You might also like

- F860 – Single Stage Compressor Test Unit
- F300 – Compressible Flow Range
- F110 – Pressure Measurement Bench

Weights & Dimensions

- 1340(L) x 900(W) x 1300(H) mm (approx.)
- Net Weight: 160 kg (approx.)
- Gross Weight: 232 kg (approx.)
- Packing Case Dimensions: 1.9(L) x 0.77(W) x 1.4(H) m
- .

All brand and/or product names are trademarks of their respective owners. Specifications and external appearance are subject to change without notice. The colour of the actual product may vary from the colour shown in the brochure.

Copyright © 2018 P.A. Hilton Limited. All rights reserved. This technical leaflet, its contents and/or layout may not be modified and/or adapted, copied in part or in whole and/or incorporated into other works without the prior written permission of P. A. Hilton Limited. Hi-Tech Education is a registered trade mark of P. A. Hilton Limited.

COUNTRY OF ORIGIN - UK WARRANTY PERIOD - 2 YEARS

Essential Services

- 380-440 Volts, Three Phase, 50Hz/60Hz, + Neutral + Earth/ground, line current up to 10A at 415v.
- 210-220 Volts, Three Phase, 60Hz + Earth/ground, line current up to 20A at 220v.
- Water: 3 litres/minute at 20m head

Ordering information

To order this product, please call PA Hilton quoting the following codes:

F865/415

F865/220

F865/415/FC

F865/220/FC

All brand and/or product names are trademarks of their respective owners. Specifications and external appearance are subject to change without notice. The colour of the actual product may vary from the colour shown in the brochure.

Copyright © 2018 P.A. Hilton Limited. All rights reserved. This technical leaflet, its contents and/or layout may not be modified and/or adapted, copied in part or in whole and/or incorporated into other works without the prior written permission of P. A. Hilton Limited. Hi-Tech Education is a registered trade mark of P. A. Hilton Limited.

COUNTRY OF ORIGIN - UK WARRANTY PERIOD - 2 YEARS